

## ITEMS SUPPLIED

Sigma 1-14 Microfuge, 24 place rotor
Microfuge tubes
Adapters (24 each)




1. Microfuge tubes of the same weight must be placed opposite each other in the rotor, or follow the scheme in 'Centrifuge Proper Balancing' diagram.
2. Adapter inserts for rotor are supplied for PCR tubes (0.2 ml tube), for 0.65 ml Microfuge tubes (use 0.5 - 0.75 ml insert). Also supplied are adapters for 0.25 - 0.4 ml tube size.
3. Ensure plastic rotor cover is in place before centrifugation.

## CENTRIFUGE Proper Balancing

The diagrams below indicate balanced loading of 2, 3 and 4 sample tubes in a centrifuge.

- Represents *sample tubes* in the holders
- Represents *empty* holders

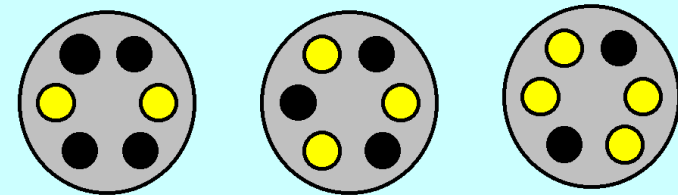

**KEEP ROTOR CLEAN  
WIPE UP ANY SPILLS IMMEDIATELY!**