

| ITEMS SUPPLIED |
|--------------------------------|
| Denver TP3002, 2 place balance |
| Weighing dishes |



1. Make sure the balance is levelled. Use the two front feet of the balance for levelling until the air bubble is centred in the level indicator.
2. Press the 'On/Off' pad to turn the balance on.
3. Place a weighing dish of appropriate size on the balance.
4. Tare the balance by pressing the 'Tare' pad.
5. **REMOVE THE WEIGHING DISH FROM THE BALANCE TO THE BENCH**

**PROPER CARE OF THE BALANCE
IS ESSENTIAL TO KEEP IT IN
GOOD WORKING ORDER**

6. Carefully transfer the material to be weighed into the weighing dish on the bench.
7. Replace weighing dish with contents on the balance and record mass.
8. Press ON/OFF pad to turn the balance off.

**CLEAN UP SPILLS
ON OR NEAR THE BALANCE
AS SOON AS THEY OCCUR**