

1. Place in freezer for at least 24 hours before use.
2. Keep the lid on when in use. The inside rows keep cooler longer than the outside rows

4°C Cooler:


This rack holds 20 microtubes (1.5 ml) and, depending on the ambient temperature, should maintain sample temperature of less than 4°C for more than 3 hours.

The rack will change from green to yellow as the temperature increases.

-20°C Cooler

This rack is a flipper rack which accommodates 20 microtubes (1.5 ml) on one side and on the opposite side 20 PCR (0.2 ml) tubes. It maintains a temperature between -20°C and -10°C for up to 3 hours

4°C COOLER

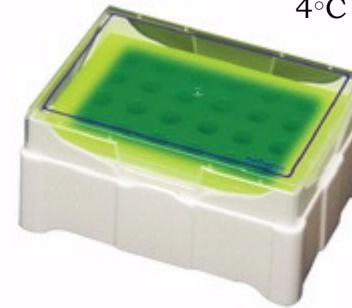

-20°C COOLER

