


UltraBasic Meter Quick Reference


1. Connect a AC adapter to the power connection.

2. Connect an electrode to the Input and ATC connectors, if desired.

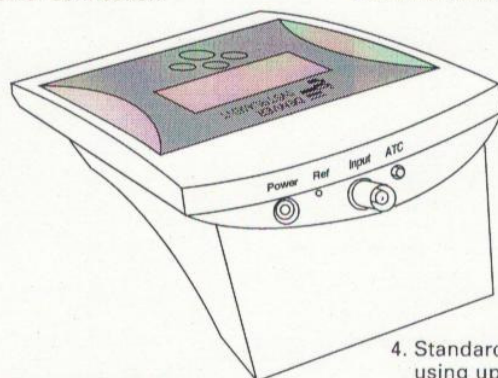

3. Press **mode** until the display indicates the appropriate measurement mode (pH or mV/relative mV).

4. Standardize the meter using up to three buffers by immersing the electrode in a buffer, stirring, then pressing **standardize** to enter each buffer.

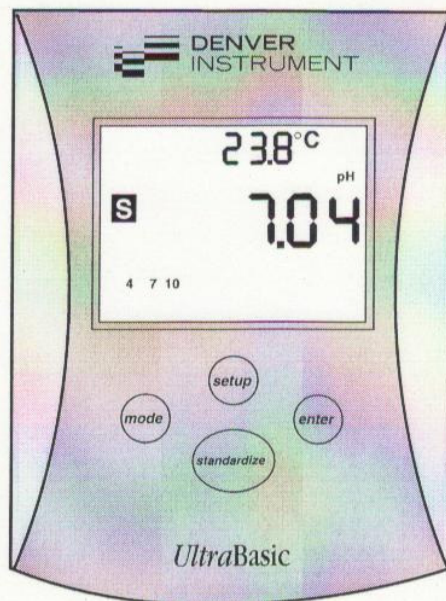

5. The display shows the current reading in pH, mV, or relative mV units.

6. Press **setup** to review electrode calibration and to clear or select buffer sets.