

<b>ITEMS SUPPLIED</b>
Benchtop Microfuge, 24 place rotor
Microfuge tubes




1. Microfuge tubes of the same weight must be placed opposite each other in the rotor, or follow the scheme in 'Centrifuge Proper Balancing' diagram.
2. Ensure plastic rotor cover is in place before centrifugation.
3. A centrifuge user manual is available at [www.kirkhoustrust.org](http://www.kirkhoustrust.org). Go to 'Resources → Research Resources → Equipment Manuals'.

## CENTRIFUGE Proper Balancing

The diagrams below indicate balanced loading of 2, 3 and 4 sample tubes in a centrifuge.

- Represents *sample tubes* in the holders
- Represents *empty* holders

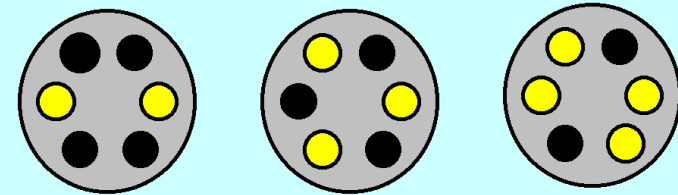

**KEEP ROTOR CLEAN**  
**WIPE UP ANY SPILLS IMMEDIATELY**